

# Annexin V-Cy5 Apoptosis Detection Kit

(Catalog #: K103-25, -100, -400; Store kit at 4°C)

## I. Introduction:

Annexin V Apoptosis Detection Kit is based on the observation that soon after initiating apoptosis, cells translocate the membrane phosphatidyl-serine (PS) from the inner face of the plasma membrane to the cell surface. Once on the cell surface, PS can be easily detected by staining with a fluorescent conjugate of Annexin V, a protein that has a high affinity for PS. The one-step staining procedure takes only 10 minutes. Detection can be analyzed by flow cytometry or by fluorescence microscopy.

## II. Kit Contents:

Components	K103-25	K103-100	K103-400
		25 assays	100 assays
Annexin V-Cy5	125 µl	500 µl	2 ml
1X Binding Buffer	12.5 ml	50 ml	2 x 100 ml

## III. Assay Protocol:

### A. Incubation of Cells with Annexin V-Cy5:

1. Induce apoptosis by desired methods.
2. Collect 1-5 x 10<sup>5</sup> cells by centrifugation.
3. Resuspend cells in 500 µl of 1X Annexin V Binding Buffer.
4. Add 5 µl of Annexin V-Cy5.
5. Incubate at room temperature for 5 min in the dark.  
Proceed to B or C below depending on method of analysis.

### B. Quantification by Flow Cytometry:

Analyze cells by flow cytometry (Ex = 649 nm; Em = 670 nm) using a Helium-Neon Laser.

For adherent cells, trypsinize and gently wash cells with serum-containing medium before incubation with Annexin V-Cy5 (A.3-5).

### C. Detection by Fluorescence Microscopy:

1. Place the cell suspension from Step A.5 on a glass slide, and cover with a glass coverslip.

For analyzing adherent cells, grow cells directly on a coverslip. Following incubation (A.5), invert coverslip on a glass slide and visualize cells. The cells can also be washed with 1X Annexin V binding Buffer and fixed in 2% formaldehyde before visualization. (Cells must be incubated with Annexin V-Cy5 before fixation because any cell membrane disruption can cause nonspecific binding of annexin V to PS on the inner surface of the cell membrane.)

2. Observe the cells under a fluorescence microscope using Cy5 filter, or FITC/Cy3/Cy5 triple band filter (Chroma Technology) if performing triple labeling using these dyes, or detect cells using a CCD camera.

Cells that have bound Annexin V-Cy5 will show bright red-blue staining on the plasma membrane.

## IV. Related Products:

### Apoptosis Detection Kits & Reagents

- Annexin V Kits & Bulk Reagents
- Caspase Assay Kits & Reagents
- Mitochondrial Apoptosis Kits & Reagents
- Nuclear Apoptosis Kits & Reagents
- Apoptosis Inducers and Set
- Apoptosis siRNA Vectors

### Cell Fractionation System

- Mitochondria/Cytosol Fractionation Kit
- Nuclear/Cytosol Fractionation Kit
- Membrane Protein Extraction Kit
- Cytosol/Particulate Rapid Separation Kit
- Mammalian Cell Extraction Kit
- FractionPREP Fractionation System

### Cell Proliferation & Senescence

- Quick Cell Proliferation Assay Kit
- Senescence Detection Kit
- High Throughput Apoptosis/Cell Viability Assay Kits
- LDH-Cytotoxicity Assay Kit
- Bioluminescence Cytotoxicity Assay Kit
- Live/Dead Cell Staining Kit

### Cell Damage & Repair

- HDAC & HAT Fluorometric & Colorimetric Assays & Drug Discovery Kits
- DNA Damage Quantification Kit
- Glutathione & Nitric Oxide Fluorometric & Colorimetric Assay Kits

### Signal Transduction

- cAMP & cGMP Assay Kits
- Akt & JNK Activity Assay Kits
- Beta-Secretase Activity Assay Kit

### Adipocyte & Lipid Transfer

- Recombinant Adiponectin, Survivin, & Leptin
- CETP & PLTP Activity Assay & Drug Discovery Kits
- Total Cholesterol Quantification Kit

### Molecular Biology & Reporter Assays

- siRNA Vectors
- Cloning Insert Quick Screening Kit
- Mitochondrial & Genomic DNA Isolation Kits
- 5 Minutes DNA Ligation Kit
- 20 Minutes Gel Staining/Destaining Kit
- β-Galactosidase Staining Kit & Luciferase Reporter Assay Kit

### Growth Factors and Cytokines

- Adiponectin/Resistin/Leptin and their Antibodies
- Recombinant Protein A and Protein G
- Recombinant Complement C5a
- Recombinant Cytokines and Growth Factors

### Monoclonal and Polyclonal Antibodies